Bedienungsanleitung User Manual

PC Software PCE-MA X

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English Contents

1	Safety notes	18
2	System requirements	18
3	Supported devices	18
4	Functions	18
5	Installation	18
6	Description of the user interface	19
7	Meaning of the individual icons on the toolbar	20
8	Operation	22
8.1	The first use of the software	22
8.2	Connect to the balance	23
8.3	Disconnect from the balance	24
8.4	Manage devices	24
8.5	Record series of measurements	26
8.6	List of measurement series	28
8.7	Numerical display	28
8.8	Statistical data	29
8.9	Graphical representation	29
8.10	Set limit values	31
8.11	Delete series of measurements	32
8.12	Export series of measurements	32
8.13	Loading and saving series of measurements	32
9	Available data	33

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1 Safety notes

Acquaint yourself with the operation of the measuring device before using it in combination with the PC software. For this purpose, use the instruction manual which comes with your PCE product. All safety notes from the device's manual have to be observed as well when the device is used with this software.

The device and the software may only be used by qualified personnel. Damage or injuries caused by non-observance of the manuals are excluded from our liability and not covered by our warranty.

We do not assume liability for printing errors or any other mistakes in this manual.

We expressly point to our general warranty terms which can be found in our general terms of business.

If you have any questions, please contact PCE Instruments. The contact details can be found at the end of this manual.

2 System requirements

- Operating system Windows 7 or higher
- USB port (2.0 or higher)
- An installed .NET framework 4.6
- A minimum resolution of 1200x600 pixels
- Optional: a printer
- Processor with 2 GHz
- 4 GB RAM recommended

3 Supported devices

- PCE-MA 50X
- PCE-MA 100
- PCE-MA 110
- PCE-MA 200
- PCE-MA 202

In chapter 9, you can find some information about how to set up the data transfer correctly as well as the data available in each case.

4 Functions

The software is used for recording, organising and evaluating series of measurements. Both weighing and drying are supported. The series of measurements can be saved to and reloaded on the PC and exported in CSV format. The data sets are displayed graphically in chronological order and also in tabular form.

The software also supports the control and monitoring of several scales. A maximum of 16 device configurations can be set up, each representing one balance. A connection can be established to all of these scales at the same time and a measurement can be carried out. Depending on the scale type, a connection via a network is also possible in addition to the connection via the COM interface.

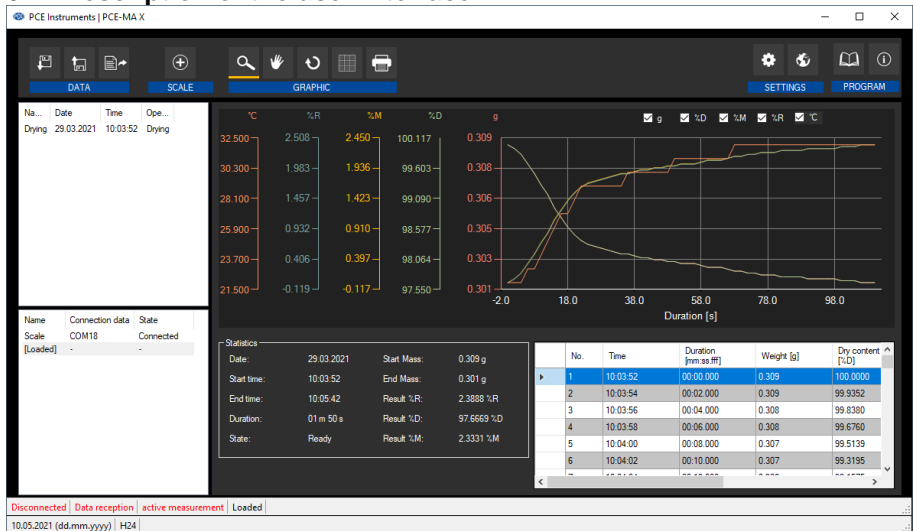
Note:

Please note that licence is required for using the software. The licence can be ordered in our online shop www.pce-instruments.com as an accessory for one of the supported devices.

5 Installation

Please run the "Setup PCE-MA X.exe" and follow the instructions of the setup.

6 Description of the user interface

The main window consists of several areas:

Below the title bar, there is a toolbar the icons of which are functionally grouped.

Below this toolbar, there is a list of measurement series (upper part) and scales (lower part) in the left part of the window.

The right part of the window is divided into an upper and a lower part.

The upper part is the area for graphical representation of the measured values.

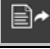
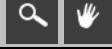
The lower part, on the other hand, provides an overview of statistical data and also a tabular listing of the measured values.

At the bottom of the main window, there are two status bars containing important information.

The lower of the two shows the static settings of the programme which can be set via a settings dialogue (date and time format and delimiter for export).

The upper status bar shows the dynamic settings of the selected balance. These include the connection status, data reception, device name, connection data, measurement settings as well as some other specific device settings.

7 Meaning of the individual icons on the toolbar

Group "Connection"	
	Connect to the selected balance
	Disconnect from the selected balance
Group "Data"	
	Create series of measurements; start recording mode
	Complete series of measurements
	Load readings from a file
	Save the readings to a file
	Export readings (CSV format)
	Delete the listed readings
Group "Balance"	
	Add new balance
	Configure selected balance
	Remove selected balance
Group "Control"	
	Request and save current reading, considering the storage conditions
	Zero balance
	Tare balance
	Set limit values
Group "Graph"	
	Enlarge the graph area ("zooming") or move the enlarged graph
	Restore the original graph
	Change the background and representation of the graph


	Print the currently visible graph
	Save the currently visible graph
Group "Settings"	
	Open settings dialogue
	Select one of the languages supported by the programme
Group "Programme"	
	Display an information dialogue
	Exit the programme

8 Operation

8.1 The first use of the software

Global software settings

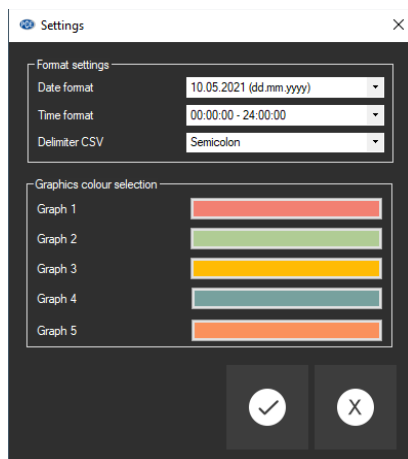Via the icon , you can open the settings dialogue. Here, you can set the date and time format. You can also select a delimiter which is used when a series of measurements is exported into a CSV file. You can also set colours for the individual graphs.

Before a balance can be connected and used with the “PCE-MA X”, it must be configured first. It is important to select the right module for the balance in the software. All modules have specific names which correspond to the product name of the balance.

Add balance

A balance can be added by clicking on the  icon.

Connection data	
Connection type:	serial interface
COM port:	COM4
Baud rate:	9600
Databits:	8
Parity:	None
Stopbits:	One

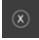
Enter the connection data of the balance in the form and save the balance configuration. Now you can connect to the balance and make all further settings for the balance and the series of measurements to be created. Further information can be found under **Configuration** 8.4.3.

8.2 Connect to the balance

When at least one balance has been configured and selected in the list of scales (see chapter 8.4.1), the connection to the device can be established by clicking on the  icon.

If the connection could be successfully established, the term "Connected" is displayed in **green** on the status bar for the dynamic data. The device settings will then be loaded and also shown in the status bar in the same colour.

8.3 Disconnect from the balance

By clicking on the  icon, an active connection to the selected balance can be terminated. If you exit the software while the connection is active, the connection to all connected devices will also be terminated.

When the software is not connected, "Disconnected" will be displayed in the status bar for the dynamic data in **red**.

8.4 Manage devices

The software allows the connection of up to 16 devices at the same time. These are listed on the left side of the software.

8.4.1 List of scales

In the left part of the main window, all scales that are currently in the software are listed.

Name	Connection data	State
Scale (0)	COM4	Disconnected
[Loaded]	-	-

Depending on the selected balance, the corresponding series of measurements are shown in the list above.

When series of measurements are loaded into the software from the memory, these are also assigned to a place marker in the list of scales.

8.4.2 Add / remove device

Click on the  button and then click on the "Apply" button to add another device to the list of scales.

To remove the configuration of a device from the software, first select a device in the list and then click on the  button to delete it.

8.4.3 Configuration

The  icon opens the configuration window of the selected moisture balance. Please note that some settings and information are only available when the software has been connected to the device.

8.4.3.1 Connection

Connection data: In order to establish a correct connection between the balance and the software, the settings of the balance (baud rate, data bits, stop bits or IP address and port) must be identical to the settings in the software.

The balance must be connected to the PC either via USB or via the network (WiFi, Ethernet).

Balance: In order for the software to communicate with the balance, the correct module must be selected. A module stands for a certain scale type.

8.4.3.2 Data

Operating mode

Depending on the type of balance, the operating modes "Drying" or "Weighing" are possible. If possible with the balance, the mode can be changed via the tab "Device settings" (see 8.4.3.3).

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To be able to call up all important information during or after a measurement, it may make sense to change the transfer method to be used as in some cases, another transfer method offers other possibilities and more information.

Two different methods of data transfer are available depending on the scale type: "Standard" and "Printer". When "Printer" is selected, data intended for a printer are transferred to the software and interpreted. This may require some extra effort when setting up the moisture balance as it is possible that not only the PC interface but also the printer interface of the device must be set.

The information on this can be called up in the software by clicking on the  icon which is located directly next to the current selection.

When a saved data record is imported, it is recommended to print via the "Printer" method. When the device is connected with the software during a live measurement, the "Standard" method is recommendable in most cases.

In chapter 9, you can find some information about how to correctly set up the data transfer with the model you have and the data available in each case. For more information, please read the manual of our device.

Storage conditions

When weighing, it is possible to save all data sets according to the sampling rate or to specify a threshold. This refers exclusively to the "main weight value". Above this threshold, the data sets are saved, otherwise they are discarded.

A tick at "Only stable values" filters out the unstable measured values.

Note:

Depending on the type of balance and method of data transfer, the software has no influence on the storage conditions so that these must be set in the balance.

8.4.3.3 Device settings

When the selected moisture balance supports further device settings, these will be listed here. By clicking on the "Apply" button, you can send the changes to the device and close the window.

8.4.3.4 Information

When the device is connected to the software, device information such as the serial number, firmware and device type will be listed.

8.4.4 Control options

The software allows you to control the moisture analyser as far as possible. You will find the most important functions in the toolbar, grouped under "Control". A click on one of the symbols sends a command to the unit.

Further functions can be found in the configuration of the respective moisture analyser under the tab "Device settings" (see chapter 8.4.3.3). The scope of the available settings differs depending on the type of device.

The following actions are possible via the buttons on the toolbar:

Request current reading, considering the storage conditions and save the reading to the active series of measurements

Zero balance

Tare balance

Set limit values. These can also be changed later for a recorded series of measurements. (see chapter 8.10)

8.5 Record series of measurements

8.5.1 Settings for the series of measurements

The settings for a measurement can be changed via the configuration dialogue of the moisture balance. These include, among others, the transfer method, the sampling rate and the storage condition. To do so, select the  icon and click on the "Data" tab. (see chapter 8.4.3.2)

8.5.2 Make a measurement

Check if the software is correctly connected to the device. To generate a series of measurements now, start recording by clicking on the  icon. In the status bar, the status of the recording is displayed.

Only then you should start weighing or drying with the balance. If the balance supports this, the data records are continuously displayed and the view is updated regularly.

During a measurement, some devices offer the possibility that the software requests data records at a set interval. This is the preferred way of measuring. Otherwise, the transfer of records is controlled via the balance:

Controlling data flow via software:

When the sampling rate has been set manually in the software, the software will request the data records from the device.

By clicking on the  button, you can additionally request the real-time value from the balance and add it to the current series of measurements. (only possible for weighing, see chapter 8.4.4)

Controlling data flow via device:

If the sampling rate option is not available in the software, the data records must be sent by the balance.

- **Print key** 

With the print key on the balance, a data set or a measurement can be sent to the software via the established connection. You may have to explicitly select "Printer" as the method for data transfer and make settings in the balance. (see chapter 8.4.3.2)
- **Continuous data transfer**

Activate continuous data transfer in your balance in order to add all data sets to the current series of measurements at the given sampling rate.

Device status:

The current status of the moisture analyzer is displayed in the status bar, e. g.

"*Connection error*" indicates a loss of connection to the meter.

"*Interpretation error*" indicates an error in interpreting the data. This may be due to incorrect software settings like e. g. an incorrect baud rate or an incorrectly selected data transfer method that does not correspond to the device. (see 8.4.3.2)

Note:

For further information regarding the possible settings for your individual device, please have a look at the data sheet of your balance.

8.5.3 Complete a measurement

To exit recording mode and complete the series of measurements, click on the  button.

8.5.4 Acoustic signals

If the balance supports the output of a signal tone sequence defined by the software, this ensures that the start of the measurement and the end of the measurement are signalled by two short consecutive tones.

8.6 List of measurement series

In the upper left part of the main window, all series of measurements belonging to the selected balance are listed. A series of measurements can be selected by a click. The tabular listing of the readings and the statistics as well as some other information will then be updated accordingly.

Name...	Date	Time	Opera...
Drying	29.03.2021	10:03:52	Drying

8.7 Numerical display

Weighing:

No.	Time	Duration [mm:ss.ff]	Weight [g]	Dry content [%D]	Water content [%M]	Wood moisture [%R]	Temperature [°C]
1	10:03:52	00:00.000	0.309	100.0000	0.0000	0.0000	22
2	10:03:54	00:02.000	0.309	99.9352	0.0648	0.0648	22
3	10:03:56	00:04.000	0.308	99.8380	0.1620	0.1623	22
4	10:03:58	00:06.000	0.308	99.6760	0.3240	0.3251	23
5	10:04:00	00:08.000	0.307	99.5139	0.4861	0.4885	23
6	10:04:02	00:10.000	0.307	99.3195	0.6805	0.6852	24
7	10:04:04	00:12.000	0.306	99.1575	0.8425	0.8497	25
8	10:04:06	00:14.000	0.306	99.0007	1.0000	1.0000	26

Drying:

No.	Date	Time	Duration [mm:ss.ff]	Weight [mg]	Net [mg]	Gross [mg]	Tara [mg]
1	29.03.2021	11:17:00	00:00.000	287.3	0.2873	287.3	0
2	29.03.2021	11:17:02	00:02.000	287.4	0.2874	287.4	0
3	29.03.2021	11:17:03	00:03.000	287.3	0.2873	287.3	0
4	29.03.2021	11:17:09	00:09.000	779.1	0.7791	779.1	0
5	29.03.2021	11:17:10	00:10.000	2347.7	2.3477	2347.7	0
6	29.03.2021	11:17:11	00:11.000	1916.4	1.9164	1916.4	0
7	29.03.2021	11:17:12	00:12.000	3586.9	3.5869	3586.9	0
8	29.03.2021	11:17:13	00:13.000	4717.7	4.7177	4717.7	0

In the lower area of the main window, all measured values of a measurement series are displayed in tabular form. Depending on the operating mode, different tabular columns are displayed.

If the balance is overloaded or if the set limit values are exceeded, the font of the relevant rows is **red**.

8.8 Statistical data

Statistics			
Date:	29.03.2021	Minimum:	279.4 mg
Start time:	11:17:00	Maximum:	12111 mg
End time:	11:17:25	Average:	5978.296 mg
Duration:	25 s	Standard devi...	4179.245 mg
Quantity:	25	Interval:	—

For weighings, the statistics window shows the time frame and some common statistical calculations.


Statistics			
Date:	29.03.2021	Start Mass:	0.309 g
Start time:	10:03:52	End Mass:	0.301 g
End time:	10:05:42	Result %R:	2.3888 %R
Duration:	01 m 50 s	Result %D:	97.6669 %D
State:	Ready	Result %M:	2.3331 %M

In drying mode, results and calculations of the currently selected drying are displayed. These include the start and end mass of the sample and the results %M (water content), %D (dry content) and %R (wood moisture).

8.9 Graphical representation


In weighing mode, the graph shows the course of the weight over time. During weighings, the limits are also marked if they have been set.

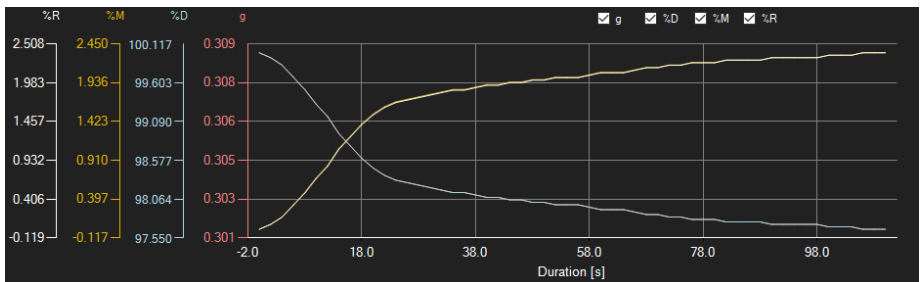In drying mode, a maximum of five measured variables can be displayed simultaneously: dry content (%D), water content (%M), wood moisture (%R) and weight (g) and temperature (°C). The tick marks at the top of the graph can be used to hide or show the corresponding graph.

Info window

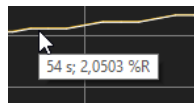Moving the mouse cursor over a dot or the curve makes a small information window with the data of the current reading pop up after a short time.

Enlargement/zooming

A freely selectable part of the displayed graph can be enlarged. For this purpose, the corresponding icon in the toolbar ("Enlarge the graph area ("zooming") or move the enlarged graph) must be a magnifying glass .

Then, a rectangle can be drawn over a part of the graph by holding the mouse button down. When the mouse button is released, the selected area appears as a new graph.

As soon as at least one enlargement has been made, it is possible to switch from enlargement mode to shift mode by clicking on the icon ("Enlarge the graph area ("zooming") or move the enlarged graph) with the magnifying glass.

Shift mode is represented by the  icon.

If the mouse is now placed over the graph area and then the left mouse button is pressed, the displayed section can be moved by holding the mouse button down.

Another click on the  icon changes back to enlargement mode, which is recognizable by the  icon.

Restoring complete view

The original graph can be restored at any time by clicking on the  icon next to the magnifying glass or hand.

Activating the background

The background of the graph and its representation can be changed via the  icon to the right. A click on the icon works like a switch: A single click makes the division of the background finer and adds some more dots to the graph. A further click on the icon changes back to standard view.

Printing the graph

The currently displayed graph can also be printed.

A click on the  icon opens the "Print" dialogue.

Saving the graph

To save the currently displayed graph, click on the  icon. The following formats are possible for the graph: BMP, JPEG, PNG, GIF, TIFF and WMF.

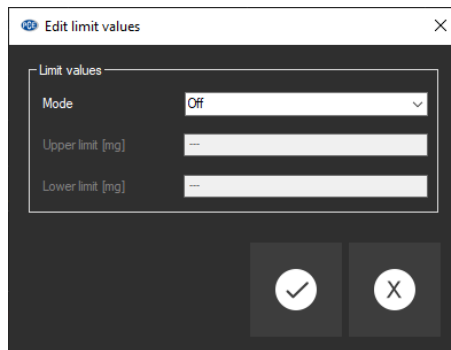
8.10 Set limit values

For each series of measurements for a weighing process, a limit mode and the corresponding limit values can be set. These will be considered when the series of measurements is saved via the  icon.

By using limit values, deviating measured values are highlighted in the chart by red, bold letters and in the graph by red horizontal lines.

To set the limits, click on the  icon in the toolbar. You can now define the modes and the corresponding limit values in the window that opens. The following modes can be selected:

1. **“Off”**
The limit values of the series of measurements are deactivated.
2. **“Upper limit”**
An upper limit can be defined. Values above the “upper limit” are outside the valid range.
3. **“Lower limit”**
A lower limit can be defined. Values below the “lower limit” are outside the valid range.
4. **“Beyond”**
An upper and lower limit can be defined. Values outside the set range are outside the valid range.
5. **“Within”**
An upper and lower limit can be defined. Values within the set range are outside the valid range.

8.11 Delete series of measurements

To delete one or more series of measurements, mark it or them in the list of measurement series and click on the  icon. Alternatively, you can use the “Del” key to do so. If there are new or modified series of measurements, the user will be asked to save these.

8.12 Export series of measurements

Via the  icon in the toolbar, the complete content of an individual series of measurements can be exported in CSV format. When the series of measurements is for a weighing process, you can also select a section from the chart (see chapter 8.7) and export it along with statistics that are adapted accordingly.

The delimiter can be changed as required via the settings dialogue .

8.13 Loading and saving series of measurements

To save a series of measurements, it must first be highlighted in the list of measurement series.

By clicking on the  button, you can now specify a file name and a target folder for this series of measurements and save it to this folder.

To load a series of measurements, you can click on the  button and then select the target folder and a file name or you can “drag & drop” the file from the Windows Explorer to the list of measurement series in the software.

Loaded series of measurements are assigned to an entry in the list of moisture balances which serves as a place marker.

9 Available data

Moisture analyser	Weighing	Drying	Configuration
PCE-MA 50X	Method "Standard"		
	Weight	Start mass End mass %M (data series) %D (data series) %R (data series) Weight (data series)	[Settings]-> [Device]-> [PC] Set up COM1, USB, Ethernet or WiFi
PCE-MA 50X	Method "Printer" [1][2]		
	Weight Net Gross Tare (Meas. settings)	Start mass End mass %M (data series) %D (data series) %R (data series) Weight (data series) Temp. (data series) (Meas. settings)	1. Additionally set up printer interface: [Settings]-> [Devices]-> [Printer] 2. Activate at least the following printout parameters: <u>Drying:</u> Start date, start time, drying parameters, start mass, drying time and result, status, separator line, end mass, end time, result <u>Weighing:</u> Weight, date, time
PCE-MA 100 PCE-MA 110 PCE-MA 200 PCE-MA 202	Method "Standard"		
	Weight (Meas. settings)	Start mass End mass %M (data series) %D (data series) %R (data series) Weight (data series) Temp. (data series) (Meas. settings)	Enable continuous data transfer (Press and hold MENU until "t" is displayed)
PCE-MA 100 PCE-MA 110 PCE-MA 200 PCE-MA 202	Method "Printer"		
	Weight Net Gross (Meas. settings)	Start mass End mass %M %D %R Weight (Meas. settings)	1) Manual printing (Press and hold MENU until "Print" is displayed) 2) Automatic printing (Press and hold MENU until "AUTO" is displayed)

^[1] The connection type (COM1, USB/ USB PC, Ethernet, WiFi) must be the same as in the configuration for the "Standard" method.

^[2] The "Printer" method is currently not available when the software is connected with the meter via the network.

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User manuals in various languages (français, italiano, español, português, nederlands, türk, polski, русский, 中文) can be found by using our product search on: www.pce-instruments.com

Specifications are subject to change without notice.

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PCE Americas Inc.
PCE Teknik Cihazları Ltd.Şti.