## PCE-PH 228

### pH-Messgerät | pH meter

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## Quick Start Guide

### Safety notes

Please read this manual carefully and completely before you use the device for the first time.

You can obtain the manual via the QR code on the front or on the supplied USB pen drive or on the respective device website at [www.pce-instruments.com](http://www.pce-instruments.com).

Damage or injuries caused by non-observance of the manual are excluded from our liability and not covered by our warranty.

- The device housing may only be opened by specialist personnel of PCE Instruments.
- Do not expose the meter to moisture.
- No technical changes must be made to the device.
- Never look into the sensor when the device is switched on, otherwise eye damage may occur.

We do not assume liability for printing errors or any other mistakes in this manual. We expressly point to our general guarantee terms which can be found in our general terms of business.

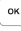
### Instructions

Press the  key to switch on the instrument.

Connect the pH electrode to the meter via the BNC connector and optionally the temperature sensor via the jack plug.

Perform a calibration before the measurement. To do this, go to the *Calibration* menu and select the buffer solutions you want to use for calibration. Then start the calibration and you will be guided through the calibration process.

Starting a calibration process is possible with the

measuring display open by a long keystroke of the  key.

Use the back button to return to the measurement screen. Now hold the electrode in the liquid to be measured. You can now read the measured value from the display.

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- Das Gerätegehäuse darf nur von Fachpersonal der PCE Deutschland GmbH geöffnet werden.