Bedienungsanleitung User Manual

PC Software PCE-PH 228

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English Contents

1	Safety notes	10
2	System requirements	10
3	Installation	10
4	Operation	11
4.1	Description of the user interface	11
5	Meaning of the individual icons on the toolbar	12
6	Operation	13
6.1	The first use of the software	13
6.2	Connect to the PCE-PH 228.....	13
6.3	Disconnect from the PCE-PH 228	14
6.4	List of measurement series	14
6.5	Numerical display	15
6.6	Statistical data.....	15
6.7	Graphical representation	16
6.8	Performing the measurement	17
6.9	Data import	18
6.10	Export of series of measurements	18
6.11	Loading and saving series of measurements.....	18

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1 Safety notes

Acquaint yourself with the operation of the measuring device before using it in combination with the PC software. For this purpose, use the instruction manual which comes with your PCE product. All safety notes from the device's manual have to be observed as well when the device is used with this software.

The device and the software may only be used by qualified personnel. Damage or injuries caused by non-observance of the manuals are excluded from our liability and not covered by our warranty.

We do not assume liability for printing errors or any other mistakes in this manual.

We expressly point to our general warranty terms which can be found in our general terms of business.

If you have any questions please contact PCE Instruments. The contact details can be found at the end of this manual.

2 System requirements


- Operating system Windows 10 or higher
- USB port (2.0 or higher)
- An installed .NET framework 4.0
- A minimum resolution of 800x600 pixels
- Optional: a printer
- Processor with 2 GHz
- 6 GB RAM recommended

3 Installation

Please run the "Setup PCE-PH 228.exe" and follow the instructions of the setup.

4 Operation

4.1 Description of the user interface

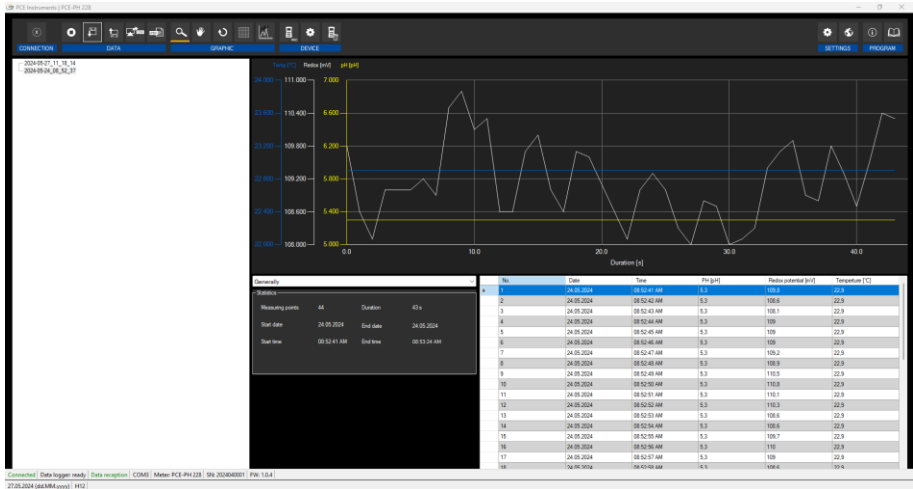The main window consists of several areas:

Below the title bar there is a "toolbar", the icons of which are functionally grouped.

Below this toolbar, there is a list of measurement series, in the left part of the window.

The right part of the window is divided into an upper and a lower part.

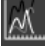
The upper part is the area for graphical representation of the measured values.

The lower part, on the other hand, provides an overview of statistical data and also a tabular listing of the measured values. At the bottom of the main window there are two "status bars" containing important information, directly above each other.

The upper status bar shows the dynamic settings of the "PCE-PH 228" which are retrieved directly from the connected device.

The lower status bar shows the static settings of the program which can be set via a settings dialog (COM port, date and time format).

5 Meaning of the individual icons on the toolbar


Group "Connection"	
	Connect to the selected device
	Disconnect from the selected device
Group "Data"	
	Create series of measurements; start recording mode
	Complete series of measurements
	Load readings from a file
	Save the readings to a file
	Export readings (CSV format)
	Create PDF report with the measured data
Group "Graph"	
	Enlarge the graph area ("zooming") or move the enlarged graph
	Restore the original graph
	Change the background and representation of the graph
	Open dialog for enabling/disabling channels and changing the channel color
Group "Device"	
	Start/Stop recording series of measurements
	Open device settings dialog
	Remove series of measurements
	Print currently visible graph
Group "Settings"	
	Open software settings dialog
	Select one of the languages supported by the programme

Group "Programme"	
	Display an information dialog
	Exit the program

6 Operation

6.1 The first use of the software

Before the software can be connected to the PCE-PH 228, the assigned COM port must be set in the software once. It can be set via the settings dialog .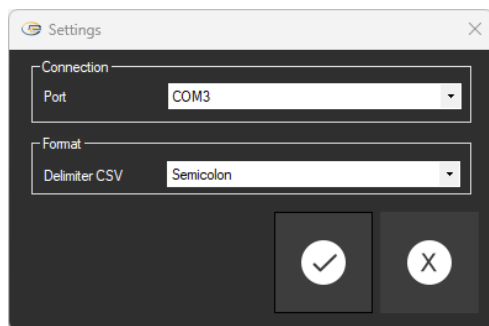

In addition to the connection data, further settings for the delimiter format for the CSV exports can be made here.

6.2 Connect to the PCE-PH 228

After the desired settings have been made and the settings dialog has been closed with a click on the "Apply" button, the connection to the PCE-PH 228 can be established by clicking on the  icon.

If the connection could be successfully established, the term "Connected" is displayed in green on the status bar for the dynamic data.

Connected	Data logger: Stopped	Data reception	COM3	Meter: PCE-PH 228	SN: 2024040001	FW: 1.0.4
27.05.2024 (dd.MM.yyyy)		H12				

If, however, no connection is established, the term "Disconnected" is displayed in red.

Disconnected	Data logger status	Data reception	COM3	Meter: PCE-PH 228	Serial number	Firmware version
27.05.2024 (dd.MM.yyyy)		H12				


6.3 Disconnect from the PCE-PH 228

By clicking on the  icon, an active connection to the PCE-PH 228 can be terminated.

If you exit the software while the connection is active, the connection will also be terminated automatically.

6.4 List of measurement series

In the left part of the main window all series of measurements, which are currently in the memory of the PC, are listed. Each series of measurements is represented in the list with the date of creation and the start time of the measurement. (e. g. 2024-05-24_08_52_37)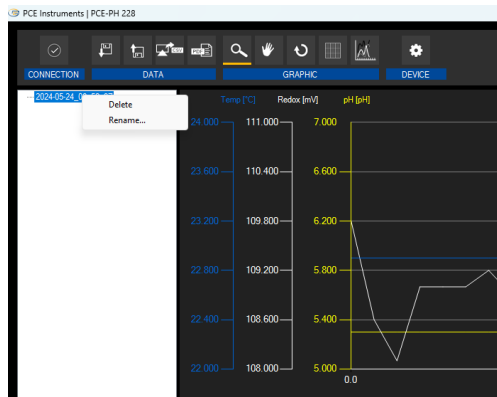

Depending on what series of measurements has been selected, the graphics, the diagram and also the statistical data are updated.

A right click allows you to delete or rename one highlighted series of measurements.

6.5 Numerical display

No.	Date	Time	PH [pH]	Redox potential [mV]	Temperature [°C]
1	24.05.2024	08:52:41	5.3	109.8	22.9
2	24.05.2024	08:52:42	5.3	108.6	22.9
3	24.05.2024	08:52:43	5.3	108.1	22.9
4	24.05.2024	08:52:44	5.3	109	22.9
5	24.05.2024	08:52:45	5.3	109	22.9
6	24.05.2024	08:52:46	5.3	109	22.9
7	24.05.2024	08:52:47	5.3	109.2	22.9
8	24.05.2024	08:52:48	5.3	108.9	22.9
9	24.05.2024	08:52:49	5.3	110.5	22.9
10	24.05.2024	08:52:50	5.3	110.8	22.9
11	24.05.2024	08:52:51	5.3	110.1	22.9
12	24.05.2024	08:52:52	5.3	110.3	22.9
13	24.05.2024	08:52:53	5.3	108.6	22.9
14	24.05.2024	08:52:54	5.3	108.6	22.9
15	24.05.2024	08:52:55	5.3	109.7	22.9
16	24.05.2024	08:52:56	5.3	110	22.9

In the lower part of the main window, all the measured values are tabulated. The chart can be sorted by columns by clicking on the respective column heading ("No.", "Time", "PH [pH]", "Redox potential [mV]" or "Temperature [°C]").

Note:

These data come directly from the PCE-PH 228; the time can thus deviate from the PC time.

6.6 Statistical data

In the general view of the statistics, you can view information about your measurement data. Here, you can see the number of measurement points, the start and end dates, as well as the start and end time of the series of measurements. You can also see the duration of measurement which is calculated from the start and end time.


General			
Statistics			
Measuring points	16	Duration	15 s
Start date	29.05.2024	End date	29.05.2024
Start time	09:56:13	End time	09:56:28

When analysing measurement data, it may be helpful to look at different views of the statistics. In addition to the general view, there are some more specific views such as of the temperature. To open this view, select the menu item "temperature" via the dropdown menu above the statistics window. Click on it to open this view.


Temperature			
Statistics			
Quantity	16	Average [°C]	22.7
Minimum [°C]	22.7	Maximum [°C]	22.7
Standard deviation	0	Total [°C]	363.2
Median [°C]	22.7	Variance [°C]	0

6.7 Graphical representation

In the area for the graphical evaluation, the selected measured values (such as Temperature, pH and Redox) are shown on the y axis and the time course (in seconds) is shown on the x axis.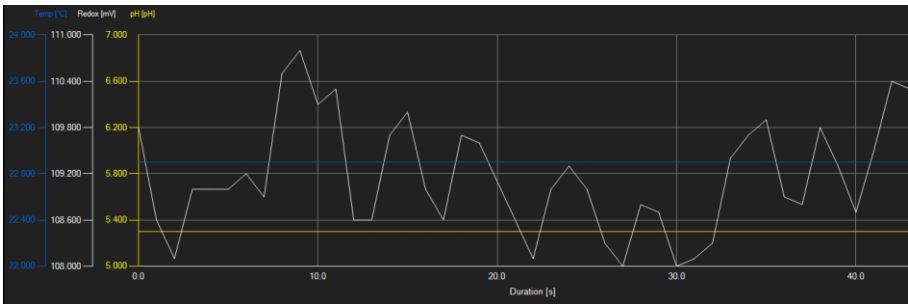

Moving the mouse cursor to a point in the displayed line makes a small information window with the data (time and unit) of the currently selected measurement value pop up after a short time.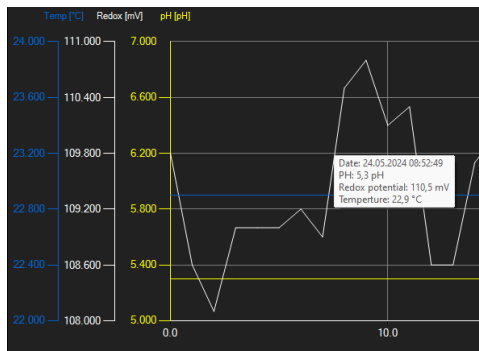

A freely selectable part of the displayed graph can be enlarged.

For this purpose, click on the  icon in the toolbar.

Then, a rectangle can be drawn over a part of the graph by holding the mouse button down. When the mouse is released, the selected area appears as a new graph.

As soon as at least one enlargement has been made, it is possible to switch from enlargement mode to shift mode by clicking the  icon.

If the mouse is now placed over the graph area and then the left mouse button is pressed, the depicted section can be moved by holding the mouse button down.


Another click on the  icon changes back to the enlargement mode.

Via the  icon to the right, some additional auxiliary lines and dots are shown in the diagram. A click on the icon works like a switch: After one click, the background is more finely divided and adds some dots to the graph. A further click on the icon changes back to standard view.

6.8 Performing the measurement

If there is an active connection to the PCE-PH 228, a new measurement can be started by clicking on the  icon.

After clicking on the corresponding icon in the toolbar ("Start a measurement"), a new dialog is then displayed, with the help of which additional options, relevant for the upcoming measurement, can be determined.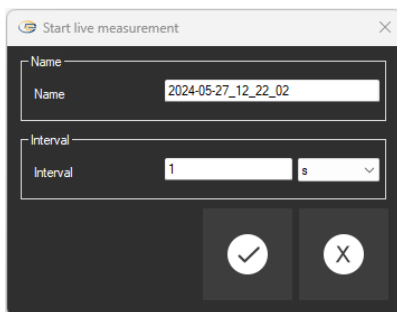


An ongoing measurement can be terminated by clicking on the  icon in the toolbar.

6.9 Data import

The internal device memory can then be read out using the software and thus the recorded series of measurements can be made available via an import.

Just click on the  icon in the toolbar. A new window will appear with the list of measurement series from which you can select which one to import.

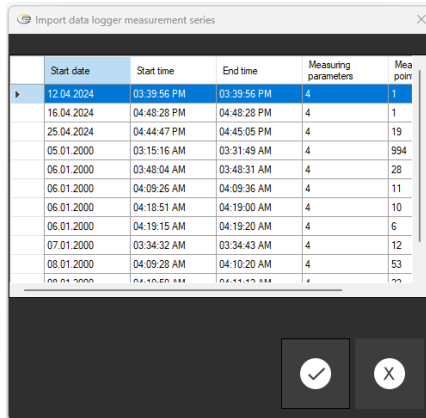This can take a few minutes, depending on the extent. The data import can now be started by clicking on the OK button or cancelled by clicking on the Cancel button.

6.10 Export of series of measurements

Measurement series can also be exported. The entire content of one measurement series can be exported in CSV format by using the  icon of the toolbar.

6.11 Loading and saving series of measurements

A series of measurements can be loaded and saved via the  and the  icon in the toolbar.

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User manuals in various languages (français, italiano, español, português, nederlands, türk, polski, русский, 中文) can be found by using our product search on: www.pce-instruments.com

Specifications are subject to change without notice.

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PCE Teknik Cihazları Ltd.Şti.