Bedienungsanleitung User Manual

PC Software PCE-SOFT SC

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English Contents

1	Safety notes	11
2	System requirements.....	11
3	Installation	11
4	Description of the user interface	12
5	Meaning of the individual icons on the toolbar.....	13
6	Operation.....	15
6.1	The first use of the software	15
6.2	Connect to the scales.....	16
6.3	Disconnect from the scales	16
6.4	List of scales	16
6.5	List of measurement series	16
6.6	Numerical display	17
6.7	Statistical data.....	17
6.8	Graphical representation	18
6.9	Control options.....	19
6.10	Performing the measurement	19
6.11	Deleting measured values	19
6.12	Exporting series of measurements	19
6.13	Loading and saving series of measurements.....	19
7	Configuration options	20
7.1	Connection.....	20
7.2	Data	20
7.3	Format	20
8	Saving data when program is closed.....	20

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1 Safety notes

Acquaint yourself with the operation of the measuring device before using it in combination with the PC software. For this purpose, use the instruction manual which comes with your PCE product. All safety notes from the device's manual have to be observed as well when the device is used with this software.

The device and the software may only be used by qualified personnel. Damage or injuries caused by non-observance of the manuals are excluded from our liability and not covered by our warranty.

We do not assume liability for printing errors or any other mistakes in this manual.

We expressly point to our general warranty terms which can be found in our general terms of business.

If you have any questions please contact PCE Instruments. The contact details can be found at the end of this manual.

2 System requirements

- Operating system Windows XP SP3 or higher
- USB port (2.0 or higher)
- An installed .NET framework 4.0
- A minimum resolution of 800x600 pixels
- Optional: a printer
- Processor with 1 GHz
- 4 GB RAM recommended


3 Installation

Please run the "Setup PCE-SOFT SC.exe" and follow the instructions of the setup.

Note:

To use the software, you need a licence (product key). The licence is related to a workplace and can only be installed on a single computer. For this software, there are two product keys which decide what functions are available: a "Basic" version (one balance) and a "Pro" version (unlimited number of balances).

4 Description of the user interface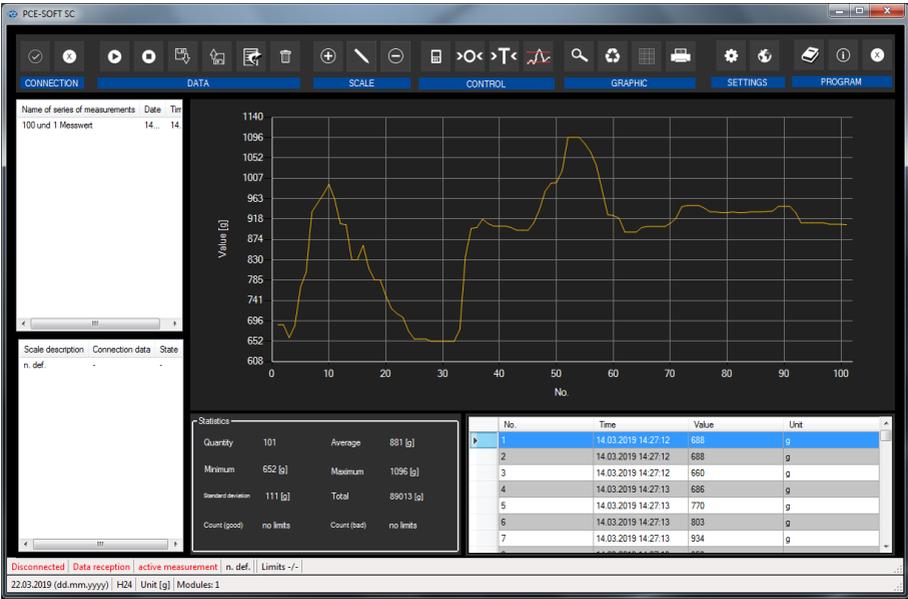

The main window consists of several areas:

Below the title bar there is a "toolbar" the icons of which are grouped according to their function.

Below this toolbar, there is a list of measurement series (above) and scales (below), in the left part of the window.

The right part of the window is divided into an upper and a lower part.

The upper part is the area for graphical representation of the measured values.

The lower part, on the other hand, provides an overview of statistical data and also a tabular listing of the measured values.

At the bottom of the main window there are two "status bars" containing important information.

The lower of the two bars shows the static settings of the program which can be set via a settings dialog (unit, date and time format).

The upper status bar shows the dynamic settings of the selected balance (COM port, set limits, data reception, selected module and the connection status to a balance).

5 Meaning of the individual icons on the toolbar

Group "Connection"		
		Connect to the selected balance
		Disconnect from the selected balance
Group "Data"		
		Load measured values/settings of a balance from a file
		Save measured values/settings of the selected balance
		Export measured values of the selected balance
		Delete measured values of the selected balance
Group "Balance"		
		Add new balance
		Configure selected balance
		Remove selected balance
Group "Control"		
		Send command to balance to get current / stable (can be selected) measured value
		Send Zero command to scales
		Send Tare command to scales
		Set limit values (will also be sent to scales, if possible)

Group "Graphics"		
		Enlarge the graphics area ("Zooming") or move the enlarged graphics
		Restore the original graphics
		Change the background and representation of the graphics
		Print the currently visible graphics
Group "Settings"		
		Open settings dialog
		Select one of the languages supported by the programme
Group "Program"		
		Display an information dialog
		Exit the program


6 Operation

6.1 The first use of the software

Before a balance can be connected to the “PCE-SOFT SC”, it must be configured. For communication of the software with the scales, it is very important to install the correct module for the scales in the software.

All modules have specific names which are related to the product names of the scales. For example, for a PCE-SD 60 CR balance, you would need the module “PCE-SD Serie”.

A balance can be added by clicking on the  icon. With the “Basic” licence, you can control one balance. With the “Pro” version, the number of scales is not limited.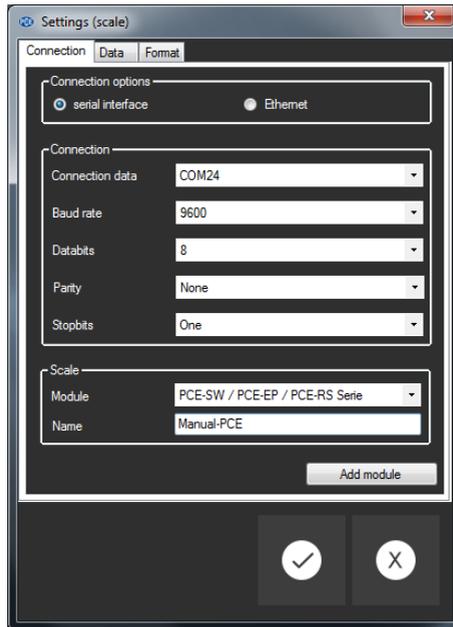

When configuring the communication with a balance, you must choose exactly the parameters appropriate for your balance.

(Serial: baud rate, data bits, stop bits, ... or Ethernet)

Further information on the configuration can be found in chapter 7.

6.2 Connect to the scales

After the desired settings have been made and the settings dialog has been closed with a click on the "Apply" button, the connection to the balance can be established by clicking on the  icon.

If the connection could be successfully established, the term "Connected" is displayed in **green** on the status bar for the dynamic data.

Connected

If, however, no connection is established, the term "Disconnected" is displayed in **red**.

Disconnected

6.3 Disconnect from the scales

By clicking on the  icon, an active connection to the selected balance can be terminated.

If you exit the software, all connections to the scales will be terminated and the configurations will be saved. (see chapter 8).

6.4 List of scales

In the lower left part of the main window, all scales currently available in the software are listed.

Scale description	Connection data	State
Manual-PCE	COM24	Disco

Depending on what balance is selected, the corresponding series of measurements will be listed above.

6.5 List of measurement series

In the upper left part of the main window all series of measurements which belong to the selected balance are listed. A series of measurements can be selected by one click. The data and statistics will then be updated, among other things (see also 6.7 Statistical data).

6.6 Numerical display

No	Time	Value	Unit
6	14.03.2019 14:27:13	903	g
7	14.03.2019 14:27:13	934	g
8	14.03.2019 14:27:13	953	g
9	14.03.2019 14:27:14	972	g
10	14.03.2019 14:27:14	994	g
11	14.03.2019 14:27:14	961	g
12	14.03.2019 14:27:14	908	g
13	14.03.2019 14:27:14	906	g
14	14.03.2019 14:27:14	830	g
15	14.03.2019 14:27:15	830	g
16	14.03.2019 14:27:15	861	g
17	14.03.2019 14:27:15	811	g
18	14.03.2019 14:27:15	786	g

In the lower part of the main window, all measured values are shown in a chart.

6.7 Statistical data


Statistics			
Quantity	101	Average	881 [g]
Minimum	652 [g]	Maximum	1096 [g]
Standard deviation	111 [g]	Total	89013 [g]
Count (good)	no limits	Count (bad)	no limits

The statistics window shows the number of measured values, the lowest and the highest measured weight, the average value, the standard deviation, the sum and – if limit values have been set – the number of values within and outside these limits.

6.8 Graphical representation

In the area for the graphical evaluation, the mass (y axis) and the consecutive number (x axis) are shown.

If limits have been set, these will also be shown (red lines).

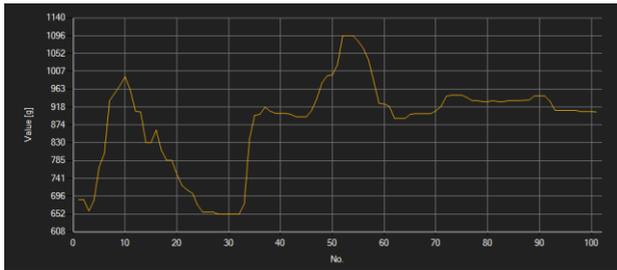Moving the mouse cursor to a point in the displayed line makes a small information window with the data of the currently selected measurement value pop up after a short time.

A freely selectable part of the displayed graphics can be enlarged.

For this purpose, the corresponding icon in the toolbar ("Enlarge the graphics area ("Zooming")

or move the enlarged graphics) must be a magnifying glass .

Then, a rectangle can be drawn over a part of the graphics by holding the mouse button down. When the mouse is released, the selected area appears as a new graphic.

As soon as at least one enlargement has been made, it is possible to switch from enlargement mode to shift mode by clicking the icon ("Enlarge the graphics area ("Zooming") or move the enlarged graphics) with the magnifying glass icon .

Shift mode is represented by the  icon.

If the mouse is now placed over the graphics area and then the left mouse button is pressed, the depicted section can be moved by holding the mouse button down.

Another click on the  icon changes back to the enlargement mode, which is recognizable by the  icon.

The original graphics can be restored at any time by clicking on the  icon next to the magnifying glass or hand.

The background of the graphics and its representation can be changed via the  icon to the right. A click on the icon works like a switch:

A single click makes the division of the background finer and adds some more dots to the graphics. A further click on the icon changes back to standard view.

The currently displayed graphics can also be printed.

A click on the  icon opens the "Print" dialog.

6.9 Control options

The control options are for easy control of the balance from the computer.

Not every balance will necessarily support all control options!

When the cursor is placed above a control element and the icon changes its colour to orange, this control element can be used.

When you click on a control element, the software will send the corresponding commands.

6.10 Performing the measurement

There are several methods to make a measurement. However, not all scales support all methods. In most cases, the settings of the balance need to be adapted first.

Information: The information about what method the selected balance supports can be found in the data sheet of the balance.

6.10.1 Keystroke on balance

Nearly any balance has a "Print" or "USB" or "RS232" key.

When this key is pressed, the balance will send the measured value.

The measured value will be received and displayed by the software (if the correct module has been selected).

6.10.2 Via software

In the "PCE-SOFT SC" software, it is possible to control the scales in order to get the current or the stable reading. To execute this command, either the  icon must be clicked or the key previously selected as shortcut key in the configuration window must be pressed first.

6.10.3 Continuously receive data from the scales

Some scales support continuous data transfer to the software if this option has been selected in their internal settings and if there is an active connection between the scales and the software.

6.11 Deleting measured values

You can always remove existing measured values. Just click on the  icon. When confirmed, the measured values/statistics will be removed and the graphics will be cleared.

6.12 Exporting series of measurements

Measurement values can also be exported. The entire content or some manually selected values from the chart can be exported in CSV format as a series of measurements by using the  icon of the toolbar.

6.13 Loading and saving series of measurements

A series of measurements and/or the PC software configuration of a balance can be loaded and saved via the  and the  icon in the toolbar.

7 Configuration options

The configuration window consists of three categories which will be explained in more detail as follows.

7.1 Connection

Connection: In order to be able to establish a correct connection between the scales and the software, the settings of the scales (baud rate, data bits, stop bits or IP address and port) must be identical with the settings in the software. The scales must be connected to the computer either via USB or via Ethernet.

Scales: To make sure the software can communicate with the scales and receive data, the correct module must be selected. The module name must be the same as the name of the scales. Some models can be used for several types of scales.

7.2 Data

Data collection: in this section, you can choose either the stable value or the current measured value. You can also select whether you want the balance to transfer data continuously. Via "raw data", you can choose that all data received from a balance are sent to output (file and/or cursor position) without checking their correctness. These functions are limited to certain scales!

Output: Here you can set how data are to be output. You can either select "Cursor position" and/or "File". If "Cursor position" has been selected, the value will be transferred from the scales to the current cursor position. If you wish to transfer the measured values to a text file, you must select the path first. How the individual fields are separated from each other is described in chapter 7.3.

Key assignment: In this area, shortcut keys can be assigned to the control elements. This function is not supported by all scales.

7.3 Format

Boxes: In this area, all important boxes can be selected. These will either be output in the file and/or at the current cursor position (Setup option see 7.2).

Delimiters: In this section, delimiters for transfer to the file and/or to the current cursor position can be set (Setup option see 7.2).

8 Saving data when program is closed

When you exit the software, the current configurations are saved to the Windows registration database and will be available again the next time the program is started.

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China

PCE (Beijing) Technology Co., Limited
1519 Room, 6 Building
Zhong Ang Times Plaza
No. 9 Mentougou Road, Tou Gou District
102300 Beijing
China
Tel: +86 (10) 8893 9660
info@pce-instruments.cn
www.pce-instruments.cn

User manuals in various languages (français, italiano, español, português, nederlands, türk, polski, русский, 中文) can be found by using our product search on: www.pce-instruments.com

Specifications are subject to change without notice.

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PCE Instruments HK Ltd.